
de	Sockel rollbar (RP)	en	Roller plinth (RP)	cs	Pojízdný sokl (RP)
el	Κυλιόμενη Βάση (RP)	en-US	Rolling stand (RP)	es	Zócalo enrollable (RP)
fr	Socle, roulant (RP)	it	Zoccolo con rotelle (RP)	nl	Verrijdbare sokkel (RP)
no	Rullesokkel (RP)	pt	Base rolante (RP)	ru	Подвижное основание (RP)
sv	Rullbar sockel (RP)				

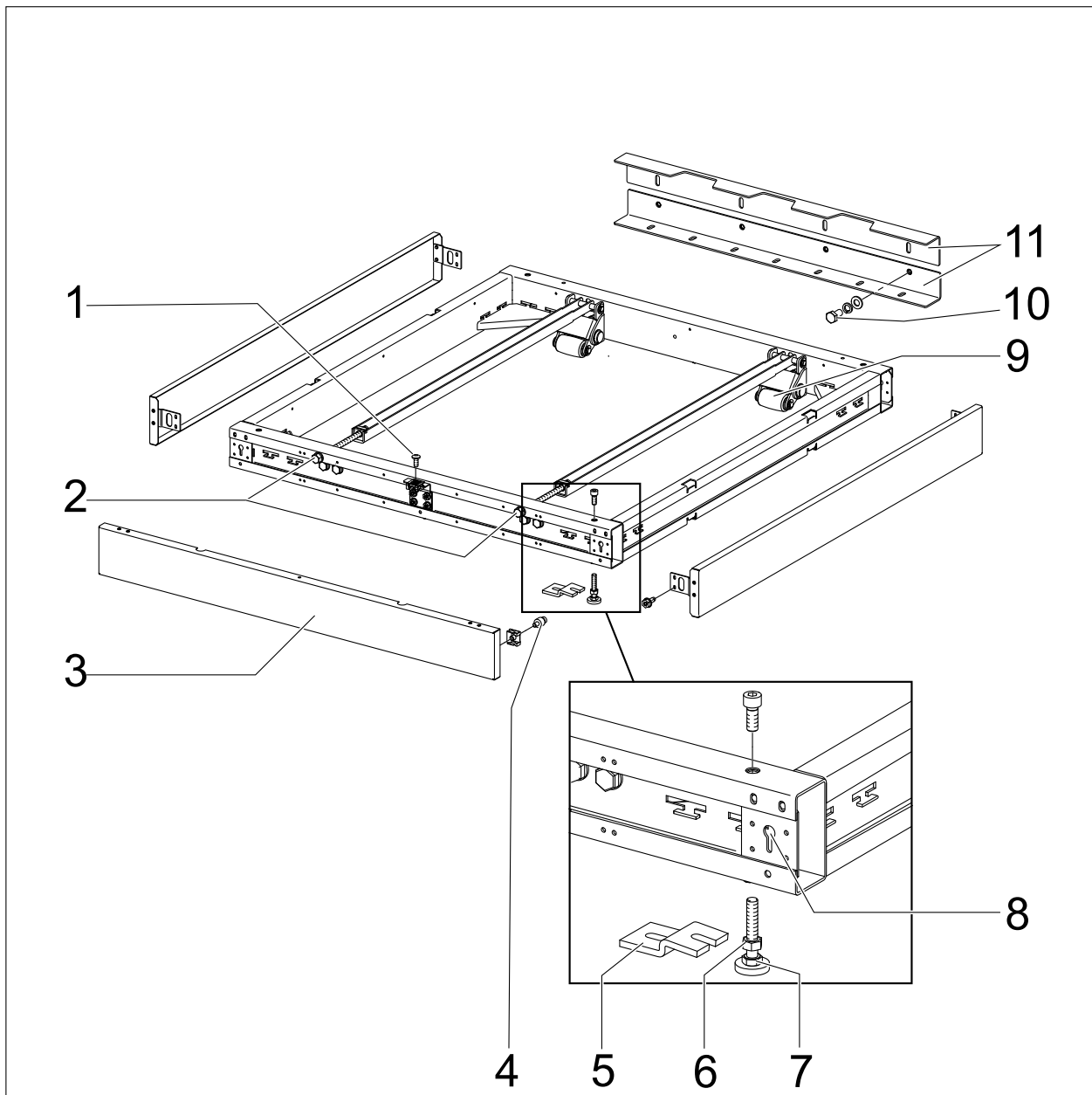1

de

Sockel rollbar (RP)

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en

Roller plinth (RP)

Reason: This conversion kit enables the roller plinth (FP) option to be fitted to PLW 8615 or PLW 8617 models.

The roller plinth (RP) enables the machine to be rolled forwards and backwards; it is not fixed to the floor.

The roller plinth (RP) can only be used with the drain pump (DP) option.

The roller plinth (RP) option cannot be combined with the automatic transport conveyor (ATC) option.

Please refer to the installation plan for the required connections, drains and infeeds.

- RP900 = suitable for 900 mm wide: PLW 8615
- RP1150 = suitable for 1150 mm wide: PLW 8617

Parts required

No.	Mat. no.	Designation
1	10941940	RP900 – Roller plinth (RP) 900 mm
1	10941950	RP1150 – Roller plinth (RP) 1150 mm

In order to carry out this work, the appropriate Technical Service Documentation (TSD) is required.

The warning and safety instructions given in the applicable model-specific TSD must be complied with.

Work must be carried out in accordance with the instructions given in the applicable model-specific TSD.

Fitting the roller plinth (RP)**Removing the plinth panel**

To level the plinth, remove the plinth panel on the infeed side (3), see Fig. 1.

- ✚ Remove the fixing screws (1) from the plinth panel (3), see Fig. 1.
- ✚ Lift the plinth panel (3) out of the keyhole brackets (8) and set it aside, taking care not to scratch it, see Fig. 1.

Levelling the plinth

The plinth has feet on the infeed side (7) and rollers on the rear (9) as well as an adjustable foot running all the way along its width (11), see Fig. 1.

- ✚ Fit the plinth and level it by adjusting the feet (7) and the adjustable foot (11) on the rear up or down, see Fig. 1.
- ✚ Once the plinth is level, tighten the counter nuts (6) on the feet (7), see Fig. 1.
- ✚ If you are installing several machines side by side, make sure that the plinth drip trays are aligned absolutely level.

Fitting optional floor mounts

The floor mounts (6) help to protect the machine in the event of an earthquake, for example, see Fig. 1.

- ✚ Mark out drill holes for the floor mount on the plinth drip tray.
- ✚ Drill holes for the plugs.
- ✚ Remove the drilling dust from the holes.
- ✚ Insert the plugs supplied with the mounts into the holes.

The special plugs supplied with the mounts will subsequently seal the drilled hole to prevent the ingress of moisture.

- ✚ Secure the floor mounts with the fixing screws.

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el**en-US****Rolling stand (RP)**

Reason: This conversion kit is for installing the optional rolling stand (RP) on models PLW 8615 and PLW 8617.

With the rolling stand (RP), the machine can be rolled back and forth; it is not secured to the floor.

The rolling stand (RP) can only be used in machines with drain pumps (DP).

The rolling stand (RP) cannot be combined with the conveyer belt option (ATC).

See the installation plan for the necessary connections, outputs and inputs.

- RP900 = suitable for 900mm-wide machines (PLW 8615)
- RP1150 = suitable for 1150mm-wide machines (PLW 8617)

Parts required

Quantity	Mat. no.	Designation
1	10941940	RP900 - 900mm rolling stand (RP = "roller plinth")
1	10941950	RP1150 - 1150mm rolling stand (RP = "roller plinth")

The TSD will be needed for all service work

The warning and safety instructions in the model-specific TSD must be observed.

Service work must be carried out following the instructions in the model-specific TSD.

Rolling stand (RP) installation**Toekick removal**

In order to align the stand, the toekick on the loading side (3) must be removed; see Fig. 1.

↻ Remove the retaining screws (1) from the toekick (3); see Fig. 1.

↻ Remove the toekick (3) from the keyhole retainers (8) and lay it down carefully to protect it from scratches; see Fig. 1.

Leveling the stand

The stand has standard feet (7) on the loading side and rollers (9) on the back side, as well as one continuous adjustable foot (11); see Fig. 1.

↻ Line the stand up and level it horizontally by tightening or loosening the standard feet (7) and the adjustable foot (11); see Fig. 1.

↻ After the stand has been horizontally leveled, tighten the counter nuts (6) on the standard feet (7); see Fig. 1.

↻ When leveling a row of machines, make sure that the drip pans are absolutely equally level.

Installing the optional floor mounting kit

The floor fasteners (6) prevent the machine from falling over, e.g., during an earthquake; see Fig. 1.

↻ Draw the drill holes for the floor fasteners on the drip pan.

↻ Drill the holes for the screw anchors.

↻ Remove any dust from the holes.

↻ Insert the provided screw anchors in the holes.

The provided special screw anchors seal the holes against any incoming moisture.

↻ Secure the floor fasteners with the retaining screws.

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